



No. 203 - Skull Bowl for the Cubic Craniophor

With damaged and fragile skulls, which cannot be properly fixed in the Cubic craniophor No. 201 with the standard skull pincers, the skull bowl is used (such as for calvaria and calotte). The skull is oriented, if at all possible, in the Frankfort Plane and fixed to the bowl with wax, plasticine, or other sticky material. The skull bowl replaces the skull pincers in the Cubic craniophor.

Caution: For cleaning please do not use acetone or acetone-containing cleaning agents.