



**No. 208 Combination instrument consisting of:**

**Mollison craniophor No. 209**

**Auricular head spanner No. 210**

The Mollison craniophor is especially suitable for adjusting the cranium in the ear-eye plane (Frankfurt horizontal).

This instrument consists of vertical metal tubes. Each bears, at its upper end, a transverse bar that can be pushed in a horizontal direction and rotated only in the axis of the vertical column. The ends of these bars are three-cornered, and the bars intended for the outer auditory canals have edges pointing upwards, whereas the edge of the third bar, which is intended to fix the lower rim of the orbit, points downwards. The instrument is collapsible as an aid to transportation.

Black's ear height gauge is fixed on to the craniophor and gives immediate readings on the millimeter scale.

**Caution: For cleaning please do not use acetone or acetone-containing cleaning agents**